



CellMosaic, Inc.
10A Roessler Road
Woburn, MA 01801, USA
Phone: 781-463-0002; Fax: 781-998-4694; E-mail: info@cellmosaic.com
www.cellmosaic.com

Thiol Assay Kit

Product Number: **CM90004**

Product Description

CellMosaic's thiol assay kit is designed to assay the free thiol group content of a biopolymer, such as a partially reduced antibody. CellMosaic routinely uses this kit for its internal bioconjugation-related research.

The assay is based on Ellman's assay using 5'5'-dithio-bis-(2-nitrobenzoic acid) (DTNB) (Ellman, G.L. 1959, Tissue sulfhydryl groups. *Arch Biochem Biophys.* 82, 70–77). Thiol is reacted with DTNB to generate 2-nitro-5-thiobenzoic acid (TNB). TNB is orange in color and has an extinction coefficient of $14,150 \text{ M}^{-1}\text{cm}^{-1}$ at 412 nm (Riddles, P.W. Blakeley, R. L., and Zerner, B. 1983, Reassessment of Ellman's reagent. *Methods Enzymol.* 91, 49-60). The amount of thiol groups will be calculated based on the amount of TNB generated.

Application of the Product

- Assay the free thiol groups.

Key Feature of the Product

- Less than 30 minutes of preparation and assay time. Fast and easy to use.

Kit Components

Four micro-centrifuge tubes per package. Each package is sufficient for 10 assays (100 μL per assay volume)

Name	Part #	Quantity
Reagent (orange label)	CM13006	1 unit
Blank (yellow label)	CM13007	1 unit
Buffer A (blue label)	CM02018	0.5 mL
Solution A (green label)	CM01003	2 mL

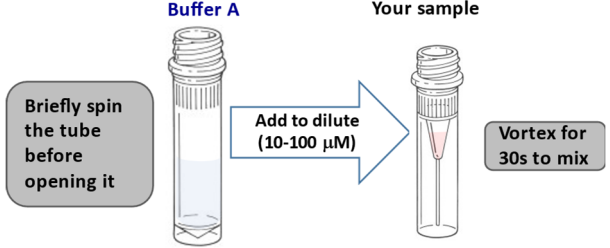
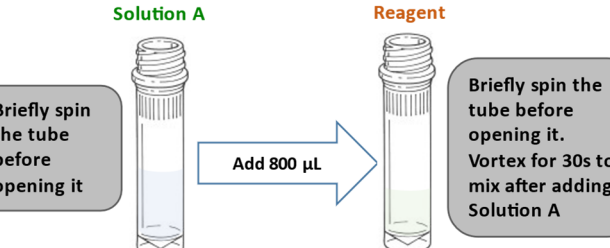
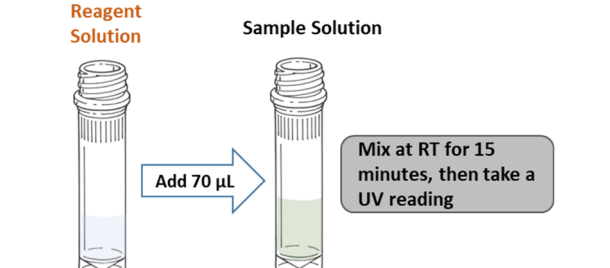
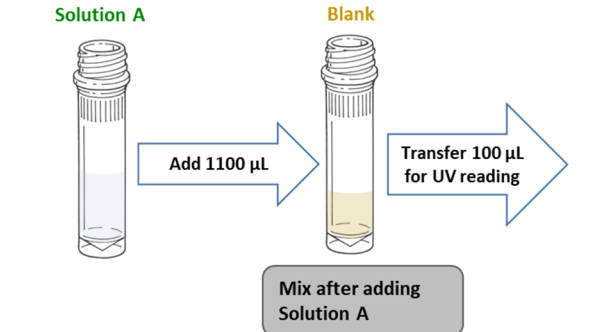
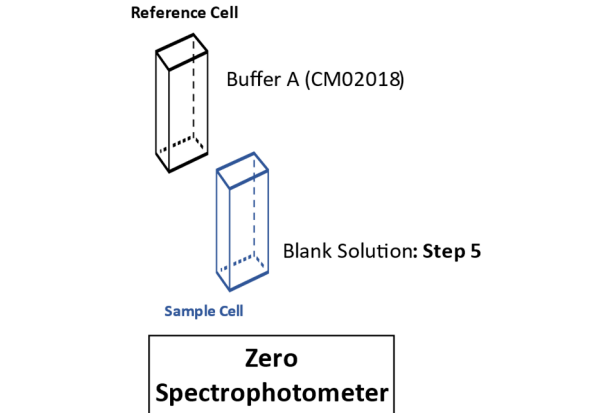
Storage/Stability

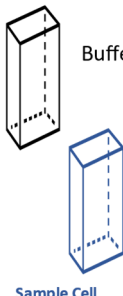
Recommended storage of the kit is at 2-8°C. For reagent and blank dissolved in solution A, they can be aliquoted and stored at -20°C up to 1 year without any sign of decomposition.

Equipment (not provided)

1. UV/vis spectrophotometer or micro-plate reader spectrophotometer with pathlength correction capability
2. Ultra-micro UV transparent cuvette: 100 μL (for UV/vis spectrophotometer) or 96-well UV microplate

Protocol

<p>1. Sample preparation: Briefly spin the tube containing Buffer A (blue label) before opening it. Dilute the biopolymer thiol in Buffer A to a total volume of 30 μL with a final concentration of thiol groups in the 10–100 μM range.</p> <p>Note: If it is an antibody with an average 4 free thiol groups per antibody, you can dilute to 1-2 mg/mL.</p>	
<p>2. Prepare Reagent solution: Briefly spin the tube containing Reagent (orange label) and Solution A (green label) before opening the tubes. Pipette 800 μL of Solution A into the Reagent tube. Vortex the solution for 30 seconds, and then centrifuge to ensure no liquid is in the cap.</p>	
<p>4. Prepare sample solution: Mix 30 μL of sample with 70 μL of Reagent solution at RT for 15 minutes.</p> <p>Note: Aliquot and store the rest of the Reagent solution at -20°C for later usage.</p>	
<p>5. Prepare Blank solution: Briefly spin the tube containing Blank (yellow label) before opening it. Pipette 1100 μL of Solution A into the Blank tube. Vortex the solution for 30 seconds, and then centrifuge to ensure no liquid is in the cap. Pipette 100 μL for the assay.</p> <p>Note: Aliquot and store the rest of the Blank solution at -20°C for later usage.</p>	
<p>6. UV readings:</p> <ul style="list-style-type: none"> • Zero the spectrophotometer with Blank solution from step 5 and Buffer A as reference at 412 nm <p>Ab (blank): ____ 0 ____</p>	

<ul style="list-style-type: none"> Measure the UV absorbance of the Sample from step 4 with Buffer A as reference at 412 nm. <p>As (sample): _____</p>	<div style="text-align: center;"> <p>Reference Cell</p>  <p>Buffer A (CM02018)</p> <p>Sample Solution: Step 4</p> <p>Sample Cell</p> <div style="border: 1px solid black; padding: 5px; display: inline-block;">Measure As</div> </div>
<p>7. Calculate the concentration of thiol</p> $\mu M = 235.5713 \times As$	$\mu M =$ _____
<p>8. Calculate the number of thiols per biopolymer (n) based on the following formula:</p> $n = \frac{\mu M (Thiol)}{\mu M (Biopolymer)}$	$n =$ _____

Important Notes & Contact Information

READ BEFORE USING ANY RESOURCES PROVIDED HEREIN

The information provided in this document and the methods included in this package are for information purposes only. CellMosaic provides no warranty of performance or suitability for the purpose described herein. Information about the chemicals and reagents used in the kit are provided as necessary.

For Research Use Only. Not for Use in Diagnostic Procedures.

The information in this document is subject to change without notice. CellMosaic assumes no responsibility for any errors that may appear in this document. In no event shall CellMosaic be liable, whether in contract, tort, warranty, or under any statute or on any other basis for special, incidental, indirect, punitive, multiple, or consequential damages in connection with or arising from this document, including but not limited to the use thereof.

NOTICE TO PURCHASER: LIMITED LICENSE

The purchase of this product includes a limited, non-transferable license to use this product to practice the labeling methods using the reagents solely for the purchaser's research activities. The license granted herein is personal to the original purchaser and may not be transferred to any other party outside the purchaser's company. No other right or license is conveyed or granted either expressly, by implication or by estoppel, to resell or repackage this product. Further information can be obtained by contacting:

Director of Licensing
c/o CellMosaic, Inc.
10-A Roessler Road, Woburn, MA 01801.
Phone: 781-463-0002
Fax: 781-998-4694
E-mail: info@cellmosaic.com